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Technical Data

- Pump head with 12 rollers (Perimax 12) provides 1 to 6 channels of liquid flow that is adjustable from 0.004 ml/min to 40 ml/min; or
- Pump head with 16 rollers phase-displaced (Perimax 16) for 1 to 3 channels of pulse-free liquid flow that is adjustable from 0.004 ml/min to 40 ml/min.
- Individually adjustable channels accommodate a wide variety of tubing materials and sizes.
- Speed range: 1 to 80 rpm, controlled by a ten-turn potentiometer with digital readout.
- High-speed key: temporarily overrides the set speed for fast-flush of the pump tubing
- Switch for off, or clockwise/anti-clockwise rotation
- Power required: 115/230 Vac at 50/60 Hz, 70 W
- Dimensions: 160x160x200 mm
- Weight: 4 to 5 kg, depending on the number of channels

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Operation

The Perimax 12 peristaltic pump is very simple to operate. The pump tubing is placed around the pump head and secured at each end in a tubing retainer. Pressure against the tubing in each channel is adjusted with a fine screw on the pressure bracket. Function keys are provided for ON/OFF and RIGHT/LEFT rotation. The rotational speed of the pump head is continuously adjustable via a 10-turn potentiometer. The digital indicator (0–999) ensures a reproducible speed setting and, therefore, a precise flow rate. A high-speed key provides for brief operation at maximum speed, for example, to obtain fast wash-out of the tubing when changing to another liquid.

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
Perimax 16 Antipuls

The 16-roller pump head consists of two half-channels of 8 rollers each in which the rollers are arranged with a 180-degree phase displacement. A piece of pump tubing is inserted into each half-channel; downstream from the pump head, a Y-fitting combines the half-channels. Consequently, a peak in one half-channel is cancelled by a trough in the parallel half-channel so that the combined flow produced is completely pulsation-free. The Perimax 16/2 can be operated with 2 pulse-free channels, 4 pulsating channels, or one pulse-free and 2 pulsating channels. Also available are the Perimax 16/1 and Perimax 16/3.

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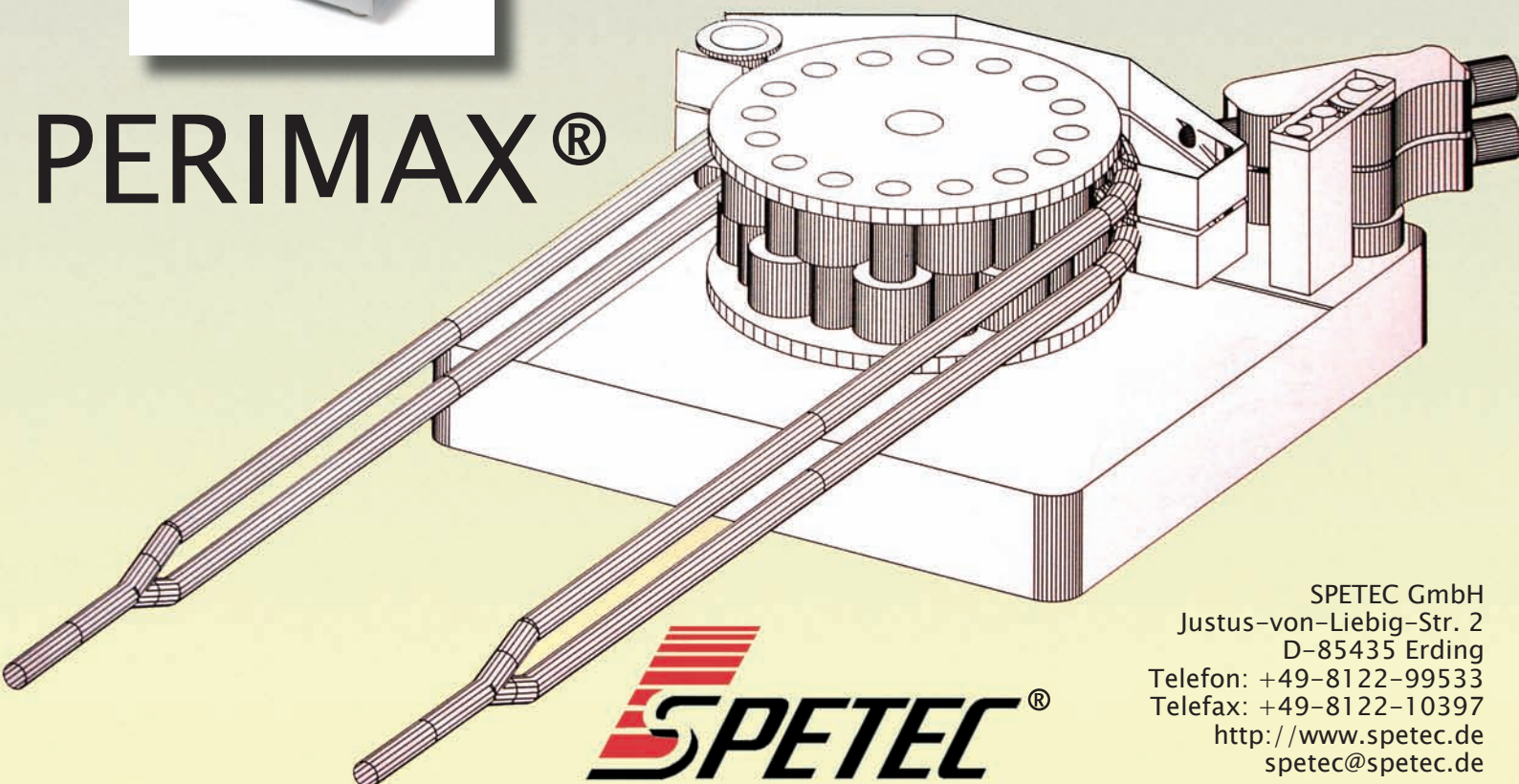Applications

The Perimax peristaltic pump is used in all branches of analytical chemistry, especially where it is essential to supply a system with accurately defined liquid quantities.

- Analytical Instrumentation
- Biotechnology, Bioanalytical Chemistry
- Liquid Chromatography
- Column Chromatography
- ICP and AA Instrumentation
- Liquid Transfer, Dosing, and Mixing
- Laboratory and Production

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PERIMAX® 16 Antipuls

drive unit

Identical drive units are employed in the Perimax 12 and Perimax 16 pumps, which differ only in the pump head. A stepper motor ensures that liquid transport is highly reproducible and constant

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Remote control

Speed and direction can be externally controlled (analog: 0-5 Vdc or digital: 26-2100 Hz) through a rear-mounted DB-9 plug.

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Tubes

Sets of pre-assembled pump tubing are available in a variety of sizes for the Perimax 16 Antipuls. A Y-fitting splits the flow between two half-channels; a second Y-fitting recombines the flows downstream.